

Anatomage

TABLE

Cadavers

How to Use the Pointer Tool

Table of Contents

How to Use the Pointer Tool

1

How to Use the Pointer Tool

Step 1: Tap the Pointer Tool icon .

Tap on a tool icon     to select the desired shape.



Step 2: Tap on the dropdown menu  to select a custom pointer color.



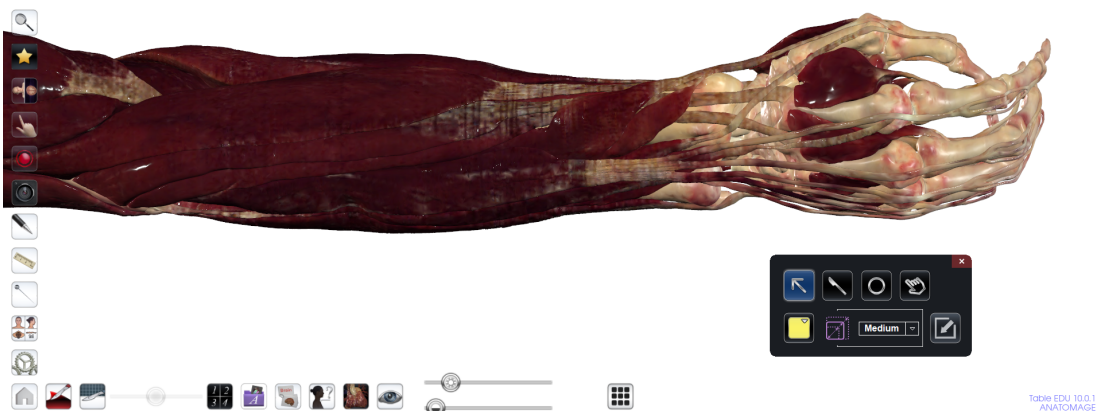


Table 10 - How to Use the Pointer Tool

Step 3: Tap on the size tab to adjust pointer tool size.

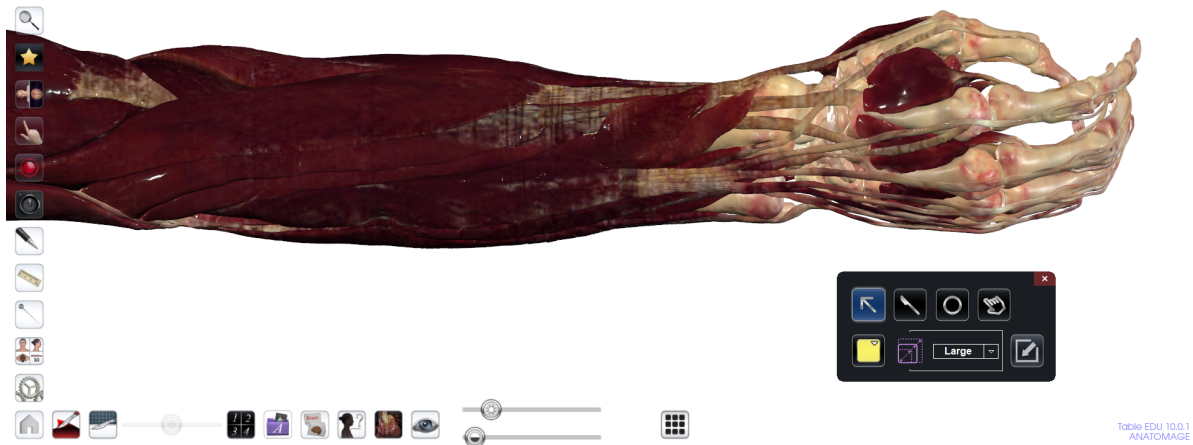

Step 4: Tap or drag your finger to move the pointer on the Table.

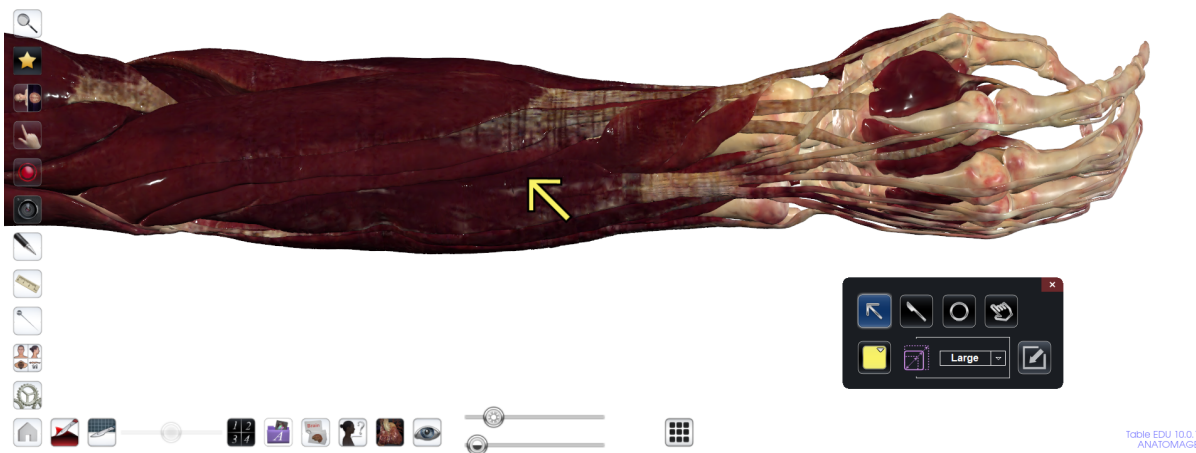



Step 5: Tap the minimize icon



to minimize the pointer dialogue. Pointer tool will

stay in use. Re-open the pointer tool by tapping the pointer icon again



.

